How to write a CD or DVD on a Linux Data Station

CD writing is no longer performed through VNMR. It is now handled by the operating system and allows the use of CDs and DVDs.

Although CD-RW and DVD-RW media can be used in principle, they entail a high risk of data loss and therefore **multi-session CD or DVD writing is strongly discouraged.** If you chose to use RW media, you do this *entirely at your own risk*. NMR staff only support CD-R, DVD-R and DVD+R and the instructions you are about to read are written for these supported media.

1. Log into your account and select the *second* desktop (Figure 1) by left clicking on the right hand side of the workspace switcher at the top of the screen (there are two desktops).

							🛛 🎾 File Browser: nmrdata		
Y				File Browser	r: nmrdata	a			-
<u>File E</u> dit <u>V</u> iew <u>G</u> o	<u>B</u> ookmarks <u>H</u> elp	_							
Back Forwar	rd Up	Stop Reload	Home Cor) 1 puter					
Location: /home1/ge	ennmr/nmrdata							🔍 50% 🔍	View as List
Information 🛩 🗙	Name	✓ Size Type	Date Modified		Owner	Group	Permissions		
~ ~	🞯 BERGENS	folder	Tue 06 Nov 2007 04	:09:43 PM MST	nmrlab	nmr	drwxr-xr-x		
	🞯 BUNDLE	folder	Tue 14 Aug 2007 10	:32:28 AM MDT	nmrlab	nmr	drwxr-xr-x		
	📁 BURIAK	folder	Wed 14 Nov 2007 0	2:26:49 PM MST	gennmr	nmr	drwxrwxr-x		
nmrdata	🎯 CAIRO	folder	Fri 17 Aug 2007 11:	02:55 AM MDT	nmrlab	nmr	drwxr-xr-x		
link to folder, 32 ite	CAMPBELL	folder	Mon 22 Oct 2007 0	:10:42 AM MDT	gennmr	nmr	drwxrwxr-x		
Sat 17 Nov 2007 09:35:1:		folder	Wed 02 May 2007 (1:10:38 AM MDT	gennmr	nmr	drwxrwxr-x		
۲	📁 ch483	folder	Thu 29 Mar 2007 01	:10:35 AM MDT	gennmr	nmr	drwxrwxr-x		
	📁 CHEM161	folder	Sat 17 Nov 2007 09	:36:35 AM MST	gennmr	nmr	drwxrwxr-x		
	📁 CHEM263	folder	Thu 18 Jan 2007 01	10:26 AM MST	gennmr	nmr	drwxrwxr-x		
	🞯 CLIVE	folder	Wed 05 Sep 2007 0	1:10:54 AM MDT	nmrlab	nmr	drwxr-xr-x		
	CORBIN	folder	Mon 29 Oct 2007 0	:49:00 AM MDT	gennmr	nmr	drwxrwxr-x		
32 items, Free space: 1	19.0 GB								
•				File Browser		Y			
32 items, Free space: 1:						<u>.</u>			
	<u>B</u> ookmarks <u>H</u> elp		Home Cor			L			
Eile Edit View Go	Bookmarks <u>H</u> elp	8 8	Home Cor	File Browser		L		§ 50% §	View as List
File Edit View Go Back Forwar	Bookmarks <u>H</u> elp	8 8		File Browser	: nmrdata		Permissions	Q 50% Q	
File Edit View Go Back Forwar Location: /home1/ge	Bookmarks Help rd Cup	Stop Reload		File Browser	: nmrdata Owner	Group	Permissions drwxr-xr-x	Q 50% Q	
File Edit View Go Back Forwar Location: /home1/ge	Bookmarks Help rd V Up ennmr/nmrdata	Stop Reload	Date Modified	File Browser	owner nmrlab	Group		Q 50% Q	
File Edit View Go Back Forwar Location: /home1/ge	Bookmarks Help The second sec	Stop Reload	Date Modified Tue 06 Nov 2007 04	File Browser puter :09:43 PM MST :32:28 AM MDT	: nmrdata Owner nmrlab nmrlab	Group nmr nmr	drwxr-xr-x	Q 50% Q	
Elle Edit View Go Back Forwar Location: /home1/ge	Bookmarks Help d V Dp ennmr/nmrdata Mame BERGENS BUNDLE	Stop Reload	Date Modified Tue 06 Nov 2007 04 Tue 14 Aug 2007 10	File Browser	: nmrdata Owner nmrlab nmrlab	Group nmr nmr nmr	drwxr-xr-x drwxr-xr-x	Q 50% Q	
Eile Edit View Go Back Fornvar Location: //home1/ge	Bookmarks Help d V Dp ennmr/nmrdata Mame BERGENS BUNDLE DURIAK	Stop Reload Size Type - folder - folder - folder - folder	Date Modified Tue 06 Nov 2007 04 Tue 14 Aug 2007 10 Wed 14 Nov 2007 0	File Browser	owner owner nmrlab nmrlab gennmr nmrlab	Group nmr nmr nmr nmr	drwxr-xr-x drwxr-xr-x drwxrwxr-x	Q 50% Q	
File Edit View Go Back Forwar Location: /home1/ge Information × × mmrdata link to folder, 32 ite	Bookmarks Help The point of the point of th	Stop Reload Size Type - folder - folder - folder - folder - folder	Date Modified Tue 06 Nov 2007 0 Tue 14 Aug 2007 10 Wed 14 Nov 2007 0 Fri 17 Aug 2007 11:	File Browser puter :09:43 PM MST :2:28 AM MDT 2:26:49 PM MST 12:25 AM MDT :10:42 AM MDT	commentations of the second se	Group nmr nmr nmr nmr nmr	drwxr-xr-x drwxr-xr-x drwxrwxr-x drwxr-xr-x	Q 50% Q	
File Edit View Go Back Forwar Location: /home1/ge Information × × mmrdata link to folder, 32 ite	Bookmarks Help The point of the point of th	 Stop Reload folder folder folder folder folder folder folder folder 	Date Modified Tue 06 Nov 2007 04 Tue 14 Aug 2007 10 Wed 14 Nov 2007 0 Fri 17 Aug 2007 11: Mon 22 Oct 2007 01	File Browser puter :09:43 PM MST :32:28 AM MDT :2:6:49 PM MST 2:55 AM MDT :10:42 AM MDT :10:38 AM MDT	Owner nmrlab nmrlab gennmr nmrlab gennmr gennmr	Group nmr nmr nmr nmr nmr nmr	drwxr-xr-x drwxr-xr-x drwxrwxr-x drwxrwxr-x drwxrwxr-x	Q 50% Q	
File Edit View Go Back Forwar Location: //home1/ge Information × × mmrdata link to folder, 32 ite Sat 17 Nov 2007 09:35:1.	Bookmarks Help The point of the point of th	 Stop Reload folder folder folder folder folder folder folder folder folder 	Date Modified Tue 06 Nov 2007 04 Tue 14 Aug 2007 10 Wed 14 Nov 2007 0 Fri 17 Aug 2007 11: Mon 22 Oct 2007 0 Wed 02 May 2007 0	File Browser puter :09:43 PM MST :32:28 AM MDT 2:26:49 PM MST 2:25 AM MDT :10:42 AM MDT 1:10:38 AM MDT	Owner nmrlab nmrlab gennmr nmrlab gennmr gennmr gennmr	Group nmr nmr nmr nmr nmr nmr	drwxr-xr-x drwxr-xr-x drwxr-xr-x drwxr-xr-x drwxr-xr-x drwxrwxr-x drwxrwxr-x	Q 50% Q	
File Edit View Go Back Forwar Location: /home1/ge Information × × mmrdata link to folder, 32 ite Sat 17 Nov 2007 09:35:1.	Bookmarks Help The provided and the pro	 Stop Reload folder 	Date Modified Tue 06 Nov 2007 0 Tue 14 Aug 2007 10 Wed 14 Nov 2007 11 Fri 17 Aug 2007 11 Mon 22 Oct 2007 0 Wed 02 May 2007 0 Thu 29 Mar 2007 01	File Browser puter :09:43 PM MST :32:28 AM MDT :2:26:49 PM MST :2:55 AM MDT :10:38 AM MDT 10:35 AM MDT 10:35 AM MST	Owner nmrlab nmrlab gennmr nmrlab gennmr gennmr gennmr	Group nmr nmr nmr nmr nmr nmr nmr nmr	drwxr-xr-x drwxr-xr-x drwxr-xr-x drwxr-xr-x drwxrwxr-x drwxrwxr-x drwxrwxr-x drwxrwxr-x	§ 50% §	
File Edit View Go Back Forwar Location: /home1/ge Information × × mmrdata link to folder, 32 ite Sat 17 Nov 2007 09:35:1.	Bookmarks Help The provided and the second	 Stop Reload - folder 	Date Modified Tue 06 Nov 2007 0 Tue 14 Aug 2007 1 Wed 14 Nov 2007 0 Fri 17 Aug 2007 11: Mon 22 Oct 2007 0 Wed 02 May 2007 0 Thu 29 Mar 2007 0 Sat 17 Nov 2007 0	File Browser puter :09:43 PM MST :32:28 AM MDT 2:26:49 PM MST 2:25 AM MDT :10:38 AM MDT 10:35 AM MDT 10:35 AM MST 10:26 AM MST	Cowner nmrlab nmrlab gennmr nmrlab gennmr gennmr gennmr gennmr gennmr	Group nmr nmr nmr nmr nmr nmr nmr nmr	drwxr-xr-x drwxr-xr-x drwxr-xr-x drwxrwxr-x drwxrwxr-x drwxrwxr-x drwxrwxr-x drwxrwxr-x drwxrwxr-x drwxrwxr-x	§ 50% §	

Figure 1: second LINUX desktop for data manipulations

2. Insert an empty CD or DVD into the DVD writer. The CD/DVD writing software, CD/DVD Creator, starts automatically and the CD/DVD Creator window appears on the desktop (Figure 2). It is best to left-click-drag this new window to the second monitor to avoid overlap with the existing windows and to make it clear which window does what. The CD/DVD Creator window should not contain any data. If data are present within the CD Creator window, a prior CD writing session may have ended ungracefully or a previous user left the check box *Reuse these files for another disc* (Figure 3) checked. In that case, select all items within the CD/DVD Creator window and move them to the

trash. Data moved to the trash from the CD/DVD Creator window will not be deleted from the hard disk or data server – only from the collection of files to be archived. But make sure that you do not drag files or directories from the two main File Browser windows into the trash. In this case data are deleted from the hard disk!

💄 🔊 🛛 🥵		🥂 🔁 🛃 🌆		File Browser: nmrdata	File Browser: CD/DVD Creator	
				📁 🎾 File Browser: nmrdata		
		File	e Browser: nmrdata			×
<u>File Edit View Go</u>						-
	, V 🔺 🧕					
Back Forwar	rd Up Sto	op Reload Home Computer				_
Location: /home1/ge	nnmr/nmrdata				🔍 50% 🔍 🛛 View as List	ž
Information 👻 🗙 🗙	Name 😽	Size Type Date Modified			CD/DVD Creator	
	DIST+FOUND	folder Sat 17 Feb 2007 01:10:31 A	File Eat View G	o <u>B</u> ookmarks <u>H</u> elp		
	DOWARY	- folder Sun 21 Oct 2007 01:10:52	м 🔍 🗸 🕨	> 🗸 🔺 🔇 🔗	6 🔊 🧭	
	MCDERMOTT	folder Wed 18 Jul 2007 09:51:08 /	Back Forw	vard Up Stop Reloa	d Home Computer Write to Disc	c
nmrdata	Mike	folder Mon 29 Oct 2007 09:49:07	Location: burn:///		Q 100	0% 🝳 🛛 View as List
link to folder, 32 ite		folder Fri 16 Nov 2007 04:43:36 P		Name 😽 Size Ty	pe Date Modified Owner Group Permis	sions
Sat 17 Nov 2007 09:35:1:	DND_083107-1-H1.fid			D_083107-1-H1.fid 4 items fol		
۲	D_083107-2-H1.fid			D_083107-2-H1.fid 4 items fol	der unknown gennmr root drwx	rwx
	D_090507-1-H1.fid			D_090507-1-H1.fid 4 items fol	er unknown gennmr root drwx	rwx
	STRYKER	folder Sun 09 Sep 2007 01:10:54		r		
	TAKATS	folder Fri 31 Aug 2007 01:10:58 A	folden Diterre			
	· · · · · · · · · · · · · · · · · · ·					
"ND_090507-1-H1.fid" s	elected		Write contents to CD			
<u>~</u>		File	Br	1 3		
<u>File E</u> dit <u>V</u> iew <u>G</u> o	<u>B</u> ookmarks <u>H</u> elp					
Back Forwar	rd Up Sto) 😵 🧑 🔛 op Reload Home Computer				
Location: /home1/ge	nnmr/nmrdata					
Information 🗸 🗙 🗙	Name 😽	Size Type Date Modified	-			
	Ø BERGENS	folder Tue 06 Nov 2007 04:09:43	PM			
	💯 BUNDLE	folder Tue 14 Aug 2007 10:32:28	АМ			
	BURIAK	folder Wed 14 Nov 2007 02:26:49	PM ^{3 items}			
nmrdata	🞯 CAIRO	folder Fri 17 Aug 2007 11:02:55 A	MMDT nmrlab nmr o	lrwxr-xr-x		
link to folder, 32 ite	📁 CAMPBELL	folder Mon 22 Oct 2007 01:10:42	AM MDT gennmr nmr o	lrwxrwxr-x		
Sat 17 Nov 2007 09:35:1.	📁 CAVELL	folder Wed 02 May 2007 01:10:38	AM MDT gennmr nmr o	Inwxnwxr-x		
1	📁 ch483	folder Thu 29 Mar 2007 01:10:35 /	AM MDT gennmr nmr o	lrwxrwxr-x		
	CHEM161	folder Sat 17 Nov 2007 09:36:35 /	AM MST gennmr nmr o	lrwxrwxr-x		
	📁 CHEM263	folder Thu 18 Jan 2007 01:10:26 A	AM MST gennmr nmr o	lrwxrwxr-x		
	💯 CLIVE	folder Wed 05 Sep 2007 01:10:54	AM MDT nmrlab nmr o	Irwxr-xr-x		
	CORBIN	folder Mon 29 Oct 2007 09:49:00	AM MDT gennmr nmr o	Inwxnwxr-x		•
	CORBIN					
32 items, Free space: 11	-					

Figure 2: CD/DVD Creator window and File Browser windows

File selection is performed in a similar manner as under MS Windows. Left click on the items that are to be removed then right click on the selected items and select *Move to Trash* from the drop down menu.

Alternatively, all items can be selected and moved to the trash by using the following menu items: Edit \rightarrow Select All Files followed by Edit \rightarrow Move to Trash.

- 3. Using the File Browser windows that are already present, navigate to the desired directory which contains data that you want to archive. Drag and drop the desired items into the CD/DVD Creator window (Figure 2). The data to be archived can be from anywhere within your nmrdata directory structure, i.e. multiple files from different directories can be archived in a single session. If an error is made, items can be safely removed from the CD/DVD Creator window by moving them to the trash (see 2, above).
- 4. Click on the button *Write contents to CD* in the CD/DVD Creator window. A new window opens as shown in Figure 3. Before writing your data, select the following:

		Write to Disc		
Write disc to : Pioneer DVD-RW DVR-112D	Information			
Disc name:	<u>W</u> rite disc to:	PIONEER DVD-RW DVR-112D		
enter a meaningful name of your choice for the CD/DVD or keep the default	Disc <u>n</u> ame:	Personal Data, Nov 20, 2007		
Data size:	Data size:	1 MB		
ensure that the data size fits on the selected media: CD: ~700 MB, DVD: ~4.7 GB	Write Options			
Write speed:	Write <u>s</u> peed:	Maximum Possible 👻		
Maximum Possible				
After Writing: ensure that the box <i>Eject disk</i> is checked and the box <i>Reuse these files for</i> another disc is <u>not</u> checked	After Writing	e files for another disc		
Click on the Write button to begin writing your CD/DVD or hit Cancel if you want to start over.	🔀 <u>H</u> elp	X <u>C</u> ancel		

Figure 3: Write to Disk window

5. Relax! your CD/DVD will be finished shortly. The writing speed depends on the media used. Do not interfere with the writing process, even though, at times, it may seem that the process is finished.

The CD/DVD will be ejected automatically once writing has completed.

 CD/DVD writing is basically a reliable process BUT it is always prudent to check whether the writing was successful.

Start VNMRJ, insert your CD, then click on the Media button (Figure 4) and navigate to the



Figure 4: The VNMRJ data manipulation buttons

cdrecorder directory by selecting it and clicking on the Set Dir button. You are now looking at the file structure on the CD/DVD. Select one *filename.fid* from the beginning of the list. Load it and process it. Repeat with a *filename.fid* in the middle and one at the end of the list. If they all process properly, you have the peace of mind that your data are on the CD/DVD in a readable format that VNMRJ can process.

If this testing sounds excessive to you, don't do it. Just remember: **the appearance of a list with filenames alone is not a guarantee by itself that the files, and not only the directory, are really on the CD/DVD**. Furthermore, these are *your* data that you want to ensure you still have when the time comes to write up your thesis or your next publication for Science or Nature!